



perform[®] sterile alcohol IPA

Sterile, alcohol based ready to use solution
for surface disinfection in cleanrooms



- Fast, effective and residue free
- Sterile (γ -irradiated) and microbial filtrated (0,2 μ m)
- Suitable for cleanrooms grade A/B, ISO 5/6
- Ready to use in a handy, sterile 500ml spray system
- Color and perfume free
- Efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to Good Manufacturing practice: Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing – current Good Manufacturing practice; U.S. Department of Health and Human Services, Food and Drug Administration, Sept. 2004

Composition

100g contain: 63,1g Propan-2-ol (equates 70% I(v/v)).

Microbiological efficacy:

Efficacy:	Concentration:	Contact time:
Bactericidal , EN1276, EN13697	Ready-to-use	1 minute
Levurocidal , EN1650, EN13697		1 minute
Fungicidal , EN1650, EN13697		15 minutes
Enveloped viruses (incl. HIV, HBV and HCV) , in accordance with (German registered association for combating viral diseases)DVV-/RKI guideline		30 sec.

Chemical-physical data:

Density	: appr. 0,88 g/cm ³ / 20°C
Color	: Colorless
pH value	: Not applicable
Flash point	: 22°C / Method: DIN 51755 Part 1
Form	: Liquid
Viscosity, dynamic	: > 2,5 mPa*s / 20°C / Raw material



Special advice

Use disinfectants safely. Always read the label and product information before use.

The sterile 500ml spray system offers the following advantages:

- Convenient size for handling and storage
- Ergonomically shaped spray pump for ease of use
- Remains sterile after opening

The bottles are sterile (with irradiation indicator) and double wrapped individually. perform® sterile alcohol IPA comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

Instructions for use

Apply perform® sterile alcohol IPA undiluted to objects and surfaces and allow to take effect. Ensure complete wetting. The product dries off quickly. perform® sterile alcohol IPA is also suitable for disinfecting gloves. Maximum application rate 50 ml/m². Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass.

Packaging unit : 10 x 500 ml bottle

Art.-no. : RREIN-PER-STIPA

hans j. michael gmbh

Gewerbegebiet Hart 11, D-71554 Weissach i.T., Phone: 0049-7191-9105-0, Fax: 0049-7191-9105-19
e-mail: office@hjm-reinraum.de, website: <http://www.cleanroom-hjm.com>